


LAQUA

Quick Manual

COND meter DS-72

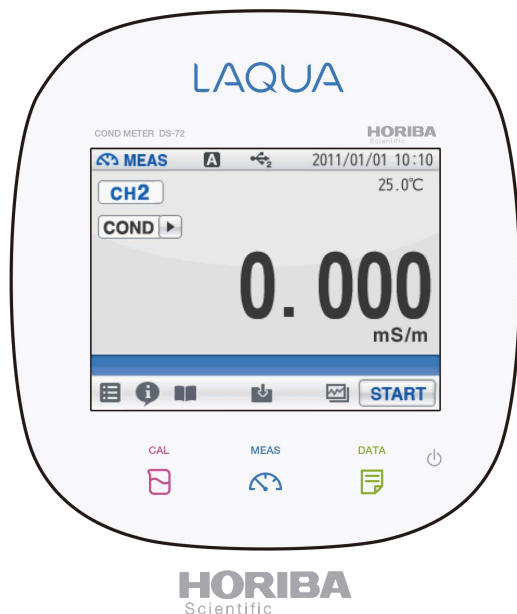


CODE : GZ0000272697A

January, 2017 © 2017 HORIBA Advanced Techno Co., Ltd.

Preparation

① Connect the electrode to the COND meter

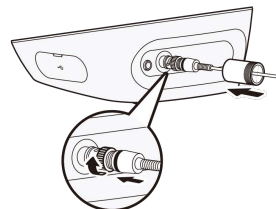

Electrode terminal



Temperature sensor terminal

② Press the power button of the COND meter




☆ Press the power button for 2 seconds.

※ Wait for 10 seconds after connecting the AC adapter before pressing the power button.

Set the cell constant

① Enter the set up mode

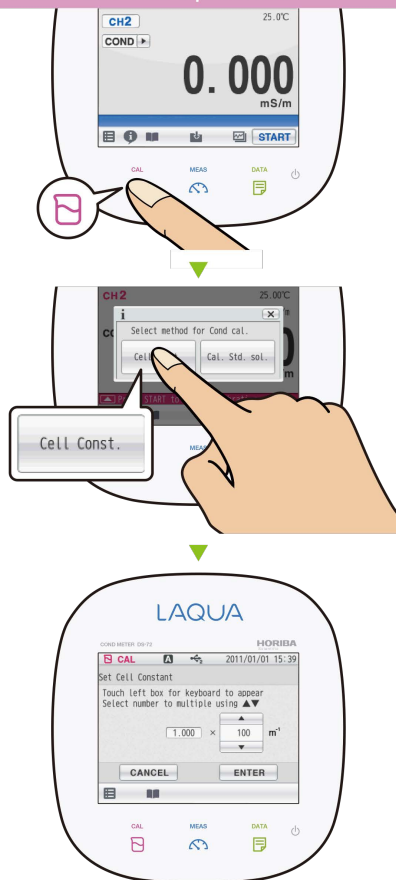


② Set the cell constant

Check the label on the electrode.

Example.)

$$3552-10D \quad 1.057 \times 100 \text{ m}^{-1} (1\text{cm}^{-1})$$

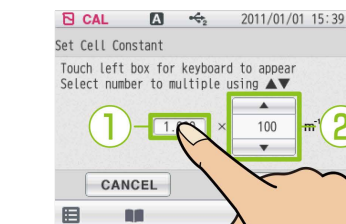

① Use the numerical keypad to enter and press the **Enter** button.




② Set the digit number with **▲ ▼** button.

Register it with **ENTER** button.

To measurement


Measurement

① Enter the measurement mode

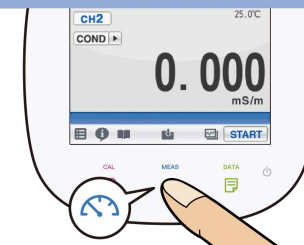


② Start the measurement

Immerse the electrode in the sample until the liquid junction part is fully submerged.

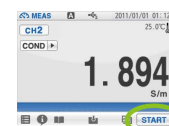

Confirm that the instantaneous values is stable, then press **START** button.

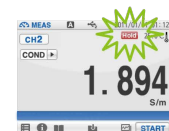

When the **Hold** indication changes from the blinking state to the lit state, measurement is complete.